

SUPACARB CP

AMAZON

HIGH PERFORMANCE CARBON BLOCK FILTER CARTRIDGE

SUPACARB CP offers the very best in activated carbon technology. The cartridge is manufactured using a unique sintered carbon that is specifically designed for maximum decolourisation and organic contaminant reduction. It offers far superior adsorption levels when compared to other carbon block cartridges.

SUPACARB CP is an extremely economical and effective solution to many common liquid treatment applications where the reduction of colour, organics and other impurities is required. The carbon block is made using a high purity micro-porous, acid washed, steam activated carbon and is particularly suited to pharmaceutical applications.

Product Features

- Unique sintered block using both granular and powdered carbon
- Spun bonded polypropylene prefiltration layer
- Modular construction from 5mm segments to 1018mm (40") long cartridges, suitable for laboratory, pilot plant and production scale operations.

The **SUPACARB CP** carbon block uses a unique sintering process that combines both granular and powdered activated carbon. This results in the highest available surface area for adsorption of contaminants. The carbon block is protected by a spun-bonded, pure polypropylene fibre prefiltration layer, to produce a unique high performance module, offering superior on-stream life.

SUPACARB CP cartridges are manufactured using the latest thermal welding technology, eliminating bypass commonly associated with adhesive bonded carbon modules. This results in a robust and durable filter assembly, particularly suited to a wide range of process conditions, including use in the harsh operating conditions encountered in pharmaceutical API plants.



SUPACARB CP FEATURES AND BENEFITS

- Combination of both granular and powdered activated carbon offers superior contaminant adsorption levels
- High active carbon content per module results in the use of smaller installations, less space required and lower costs
- Convenient low-weight cartridge format is easy to handle and modules can be changed-out without using lifting gear, special tools or other apparatus
- High purity micro-porous steam activated powdered carbon is particularly suited to removal of coloured impurities from pharmaceutical product and solvent streams.
- Graded density prefiltration layer protects carbon block offering extended life
- Thermally bonded construction eliminates fluidising, channelling and bypass
- All polymeric support structure offers excellent chemical compatibility
- Each cartridge is embossed with an identification number and fully traceable

SUPACARB CP

TECHNICAL DATA

Dimensions

Outside Diameter: 66.5mm
Inside Diameter: 27mm
Length: See ordering guide

Sterilisation and Sanitisation*1

Steam: 121°C for 15 mins (20 cycles)
Hot Water: 90°C for 30 mins (0.2 bar Δp max)

*1Applies to single open end cartridges only. For all steaming and hot water applications, the stainless steel end cap insert or Glass Filled end cap option must be used .

Maximum Operating Conditions

Temperature: 80°C
Recommended Maximum Differential Pressure: 1.5 Bar

Product validation guide available on request. All **SUPACARB CP** cartridges are manufactured under strict control with batch number identification, giving full traceability on all components.

INDUSTRIES AND APPLICATIONS

- | | |
|-----------------------------------|---|
| Pharmaceutical and Fine Chemicals | • Removal of coloured impurities from product and solvent streams |
| Food and Beverage | • Colour and odour reduction in beverages |
| Water Treatment | • Removal of chlorine, ozone, odour, colour and trace organic compounds |
| Chemical Manufacturing | • Removal of catalysts and wastewater treatment |

AMAZON FILTERS LTD.

ALBANY PARK ESTATE, FRIMLEY ROAD,
CAMBERLEY, SURREY, GU16 7PG
ENGLAND

Tel: +44 (0) 1276 670 600

Fax: +44 (0) 1276 670 101

Email: sales@amazonfilters.co.uk

Web: www.amazonfilters.co.uk



Certificate No. FM39584

Distributed by: